## REGLO-E Digital

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## Important information

### ▶ Safety

The installation of this panel-mounted pump must be carried out by a skilled person who is aware of the potential hazard involved in handling electrical components.

**Caution:** rotating pumphead creates a pinch and crush hazard.

### ▶ Wiring

See figure on front page as well as the attached wiring diagram.

### ▶ Active cooling system

Depending on the installation and the operating time, an active cooling system (ventilator) must be installed.

### ▶ Connecting cable

- Shielded cables are to be used between control unit and drive. The shield must be connected to ground on both ends.  
- The maximum length of the cable checked and approved by Ismatec is 2 meters. Depending on the situation, longer cables can be used, which however may cause malfunction. For more information, please do not hesitate to contact us.

### ▶ CE conformity

- Our standard pump models comply with the applicable EC directives.  
- However, within the EC, our panel-mounted pump models must not be operated before the complete system or instrument complies with all the corresponding EC directives.

## Technical documentation

Please refer to the enclosed:

- ▶ dimensioned diagrams
- ▶ wiring diagrams
- ▶ operating instructions for:
  - fuse ratings and voltage
  - DIP-switch settings
  - how to operate the pump

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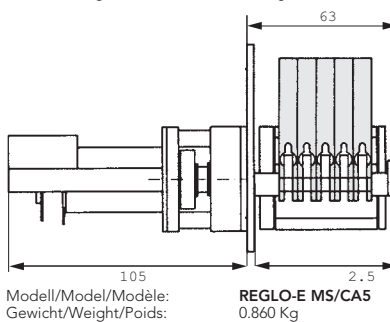
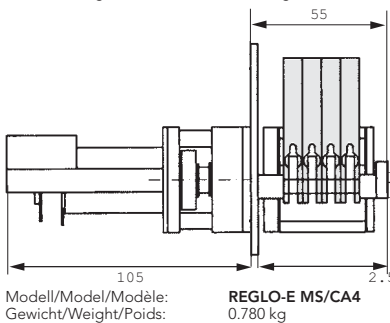
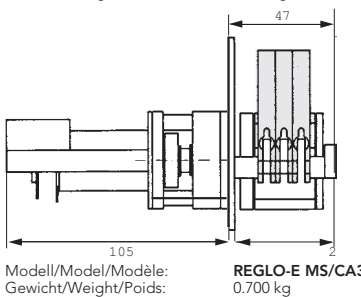
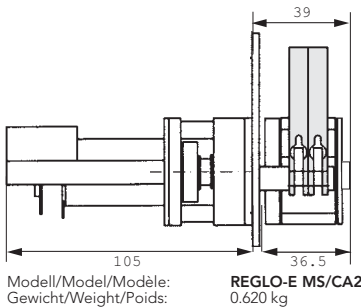
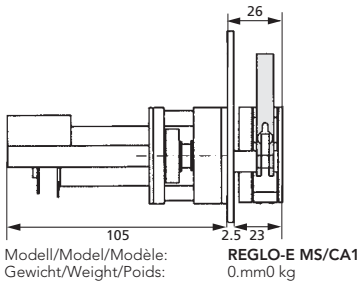
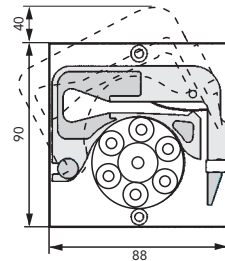
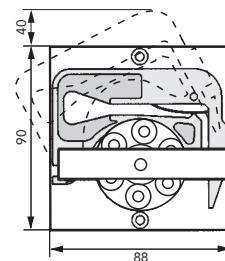
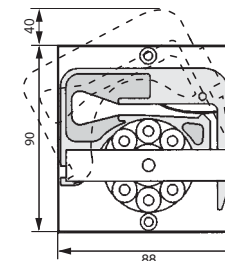
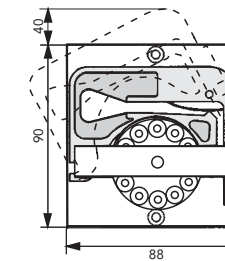
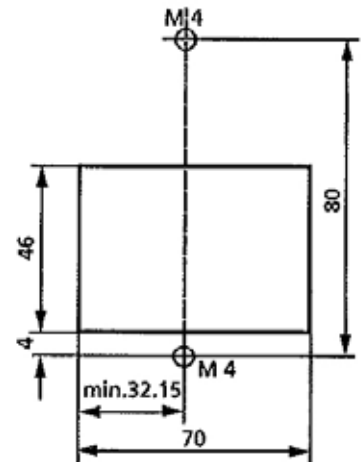
## Type of cassettes

### MS/CA Click'n'go

Spring-loaded cassette with stainless steel pressure spring providing defined and reproducible occlusion conditions. Independent of tubing diameter and material, the occlusion is automatically correct.

### MS/CA Pressure lever

Cassette with pressure lever for adjustable occlusion (may be an advantage for elevated differential pressure conditions).

## Cassette material

### POM-C

- (Polyoxymethylene-Copolymer) (Polyacetal, Polyformaldehyde)
- Good chemical resistance to many organic solvents and strong alkaline chemicals.
- Is affected by strong acids and oxidizing substances.
- UV-stabilized and stable at (dry) temperatures up to 80°C (in continuous use) or 136°C (for a short time).

### PVDF<sup>1)</sup>

- (Polyvinylidene fluoride)
- Very good chemical resistance to acids and most aliphatic, aromatic and chlorinated solvents.
- Not suitable for long contact with esters, ketones, amines and strong alkaline chemicals.
- Stable to UV and temperatures up to 110°C (continuous use) or 142°C (for a short time)

<sup>1)</sup> only cassettes with pressure lever

## Pump tubing

These cassette pumps take tubing with 3 colour-coded collars. Thanks to these collars the tubes are always inserted under the same conditions, which results in a high repeatability of the flow rate.

What is more, the different colour codes help to easily recognize the various tubing diameters.

Tubing with 3 collars can be used twice. Thanks to this advantage the service life of the tubes can be doubled.

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
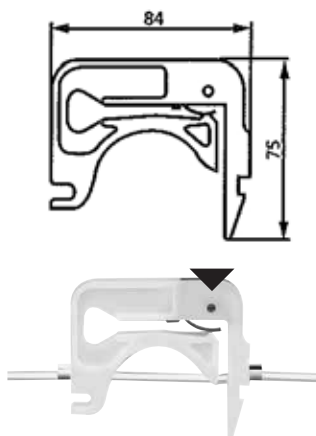
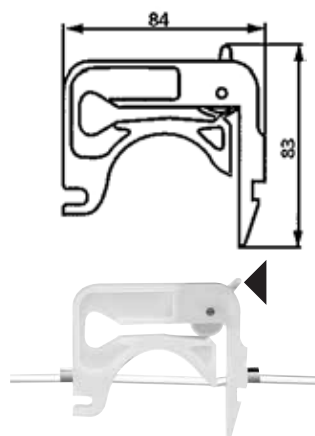
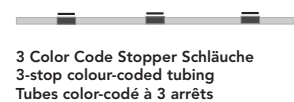
**MS/CA Anpresshebel**  
**MS/CA pressure lever**  
**MS/CA levier de pression**

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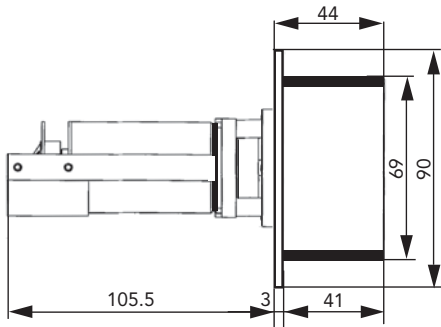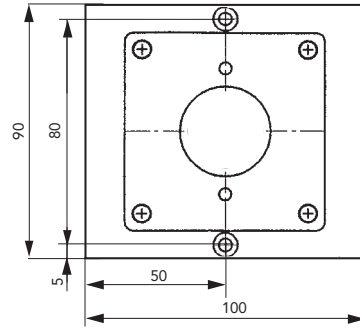
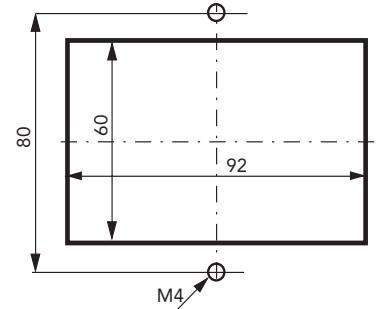
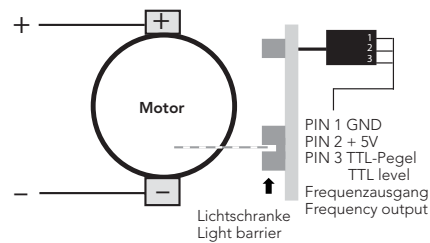
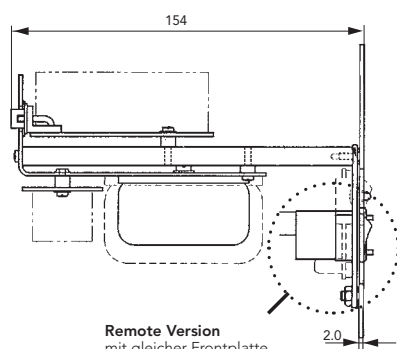
**Remote version**  
with same front panel  
but without control panel

## Installation and electrical connections

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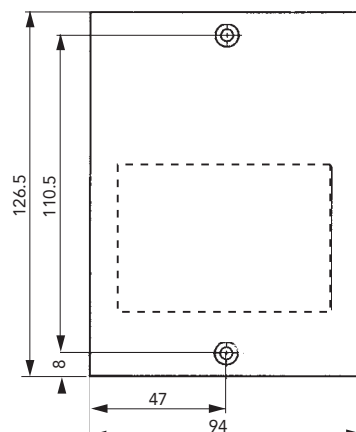
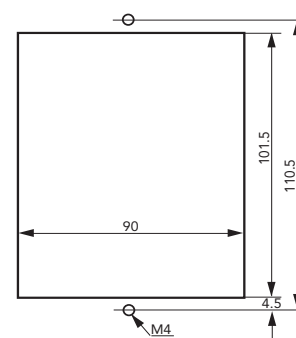
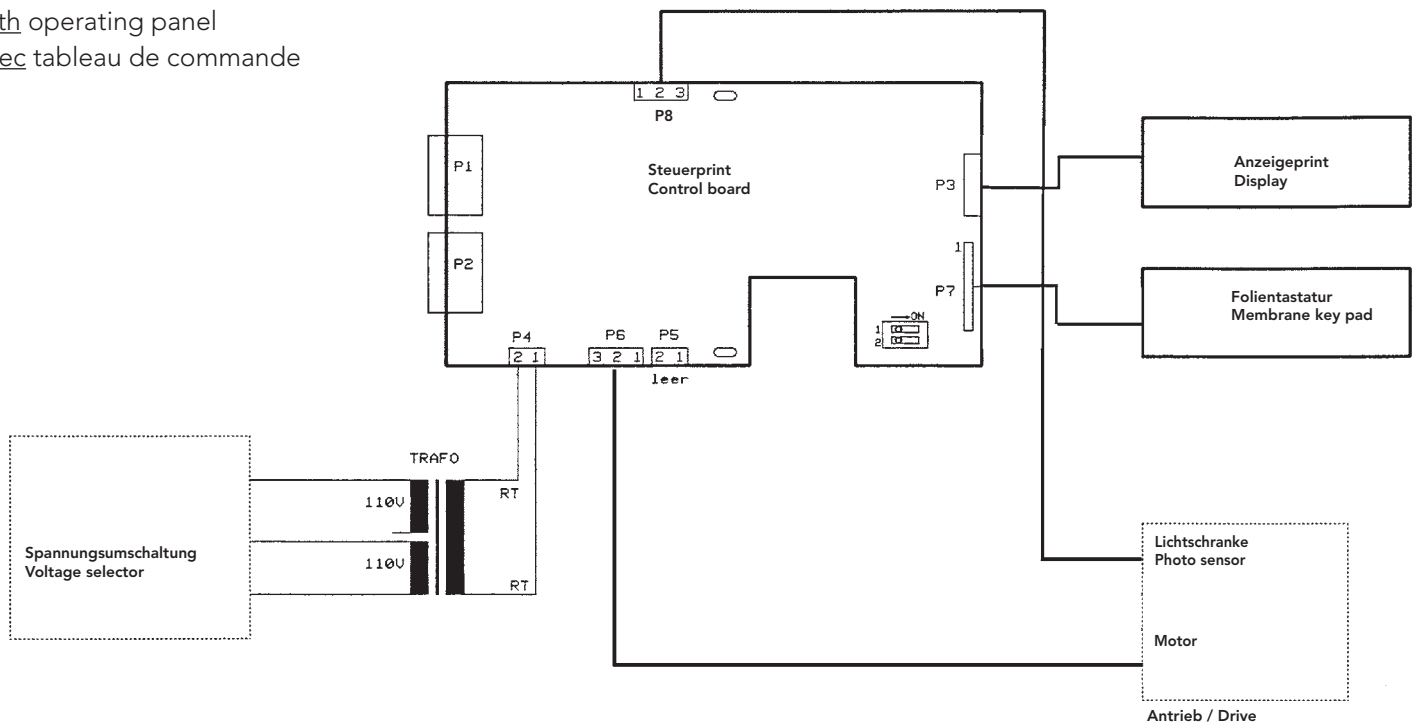
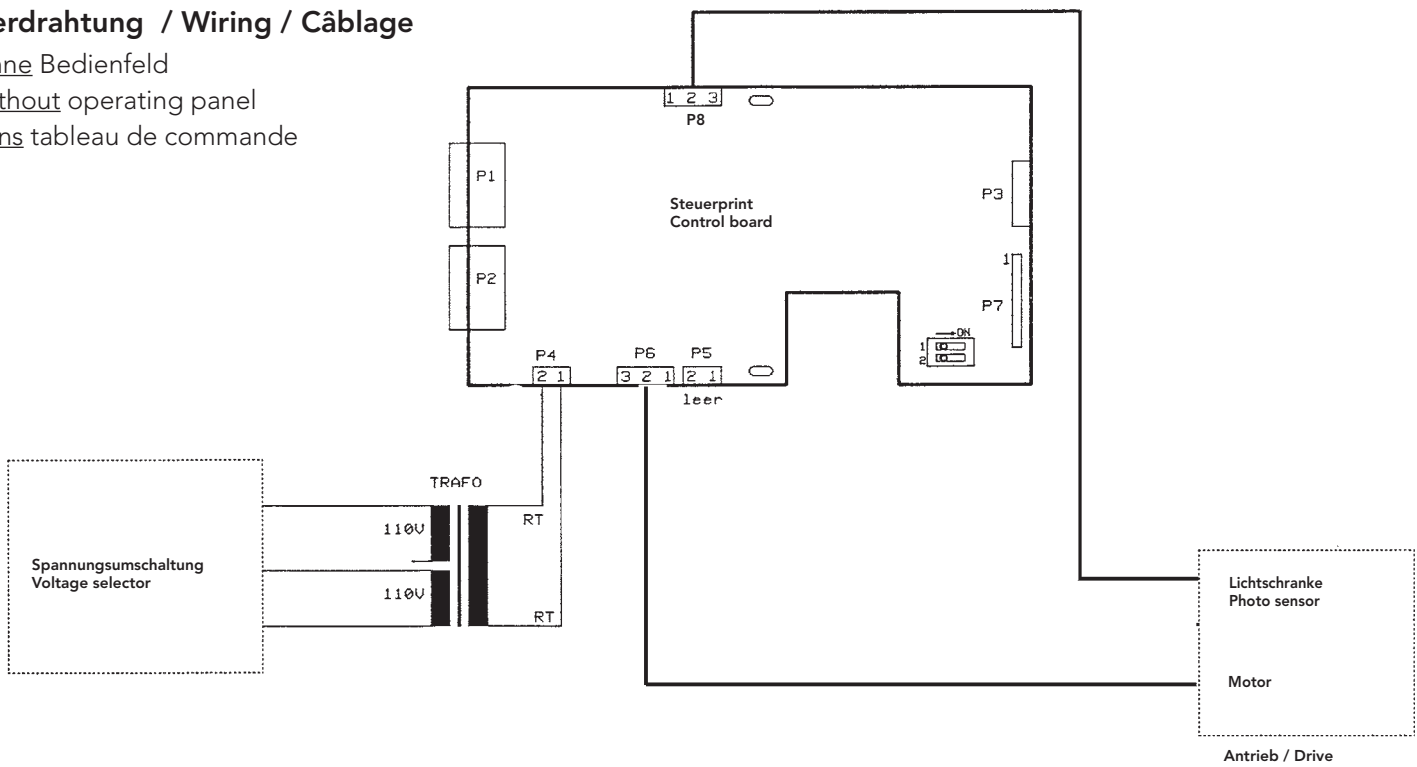
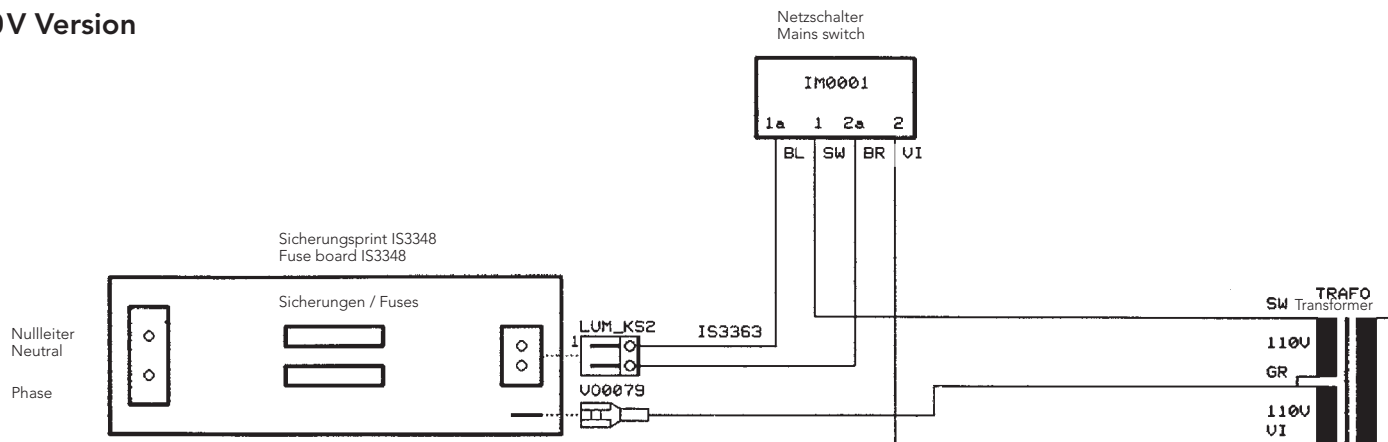
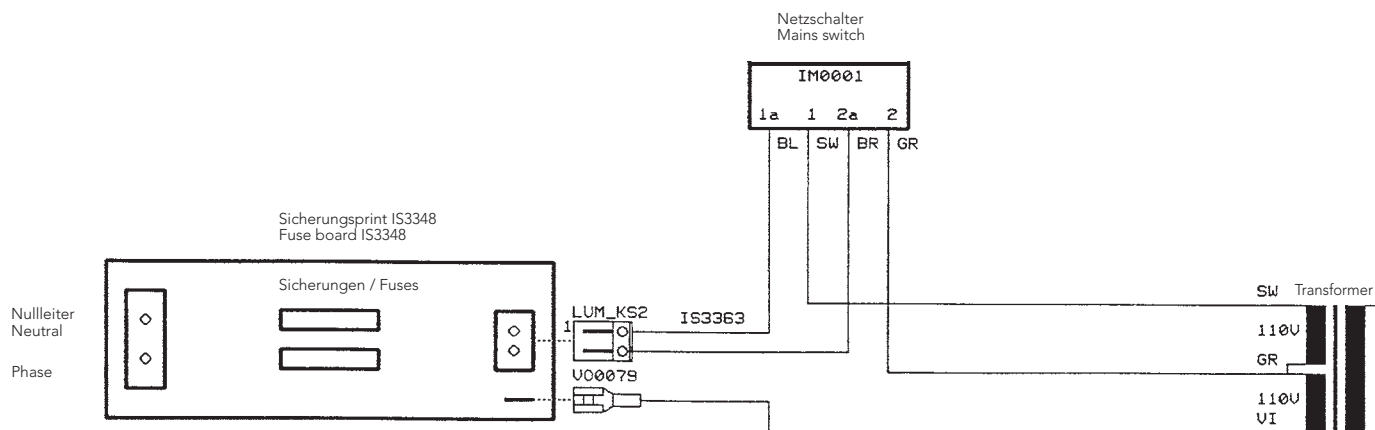
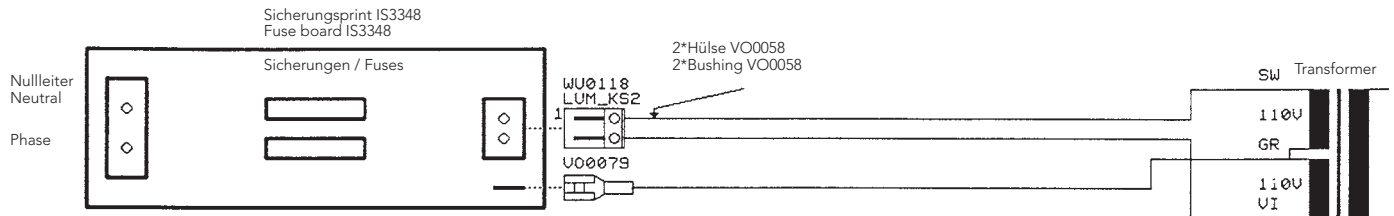
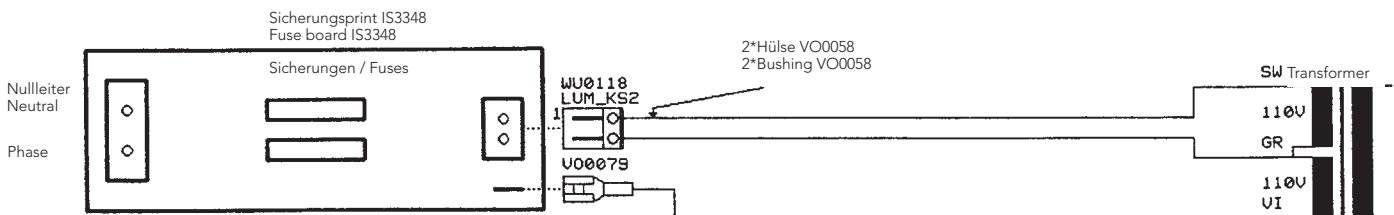
115V Remote Version

**ISMATEC®**

Learn more  
about the  
Ismatec® product  
line by visiting:  
[www.ismatec.com](http://www.ismatec.com)

**For ordering and technical support, please contact:**

**North America** [sales@ismatec.com](mailto:sales@ismatec.com) | 1-800-323-4340 | 1-847-549-7600

**Europe** [sales.europe@ismatec.com](mailto:sales.europe@ismatec.com) | +49 (0) 9377 9203-0

 **Cole-Parmer®**